

Ε

aliquot) by mixing TEV protease into protein solution (pipet well to mix thoroughly) and leaving the reaction at 4 °C undisturbed (still), WITHOUT rocking, for ~16 hours or overnight. Of course, it would be recommended to perform a test of different time-points, as it may depend from case to case, depending on many things, including TEV site accessibility.

- 1. Waugh, D.S., Protein Expr. Purif., **80(2)**, 283-293 (2011).
- 2. Cesaratto, F. et al., K. Biotechnol., **231**, 239-249 (2016).
- 3. Nallamsetty, S. et al., Protein Expr. Purif., **38(1)**, 108-115 (2004).
- 4. Zdanov, A.S. et al., Russ. J. Bioorg. Chem., **29(5)**, 415-418 (2003).